

Biocontam-Zero

Disinfection Solution for Water Bath in CO₂ Incubators

As an effective antimicrobial, SolMate™ Biocontam-Zero prevents contamination of sterile water within CO₂ incubator water baths by mycoplasma, bacteria, viruses, fungi, and yeast while preserving water quality.

Excellent antimicrobial performance

- Prevention of biological contamination through step-by-step antimicrobial activity.
- Sterilized effects for 2~4 weeks.

Non-toxic nature

- Based on sodium benzoate (Sodium benzoate, C₇H₅NaO₂), a natural compound found in certain fruits and spices.
- Non-toxic, non-corrosive, and non-volatile.



Comprehensive antimicrobial spectrum

- Targeting mycoplasma, bacteria, viruses, fungi, yeast and more

How to Use

- Allow to thaw to room temperature prior to use.
- Gently swirl the solution in the bottle (Wipe the outside of the bottle with a disinfectant solution, such as 70% ethanol, before opening)
- Recommended dilution : 1:100 in sterile water.

Ordering Information

Brand	Product	Cat. No.	Size	Conc.	Storage
SolMate™	Biocontam-Zero	SM-B01-010	100 ml	100 X	-20°C

Memo

Biocleaner-S

for Disinfection Surfaces

The ultimate solution for preventing biological contamination on incubators, laboratory equipment, and work surfaces. Trust SolMate™ Biocleaner-S to keep your lab surfaces pristine and protected from mycoplasma, bacteria, viruses, yeast, mold, and more.

Strong Anti-bacterial Effect

- Superiority over ethanol and other common lab alcohols

Easy to Use

- Convenient spray form for ease of use and storage.
- Room temperature storage

Non-toxic Materials

- Harmless chemicals to the human body
- Non-toxic, biodegradable, and non-corrosive.

Comprehensive antimicrobial spectrum

- Targeting mycoplasma, bacteria, viruses, fungi, yeast and more



How to Use

1. Apply an ample amount of Biocleaner-S. to thoroughly saturate the designated areas for disinfection, including workstations, laboratory equipment, and incubators.
2. Allow the sprayed solution to air-dry naturally without wiping it off.
3. Spray incubators once every 2 weeks.

Ordering Information

Brand	Product	Cat. No.	Size	Storage
SolMate™	Biocleaner-S	SM-B02-100	1000 ml	Room temperature